

What to do in case of eye splash

Manual for Eye washing

Diphoterine eyewash is now installed throughout the Genopode

Where?

- In each lab in the corner of the building, on Floors 3,4 and 5 near the laboratory sink a flask of Diphoterine is installed.

These labs are indicated with the sign



- In floors 1 and 2: in laboratories and premises in which chemicals are used.

Why ?

In case of chemical splash in the eyes.

Diphoterine is a solution that **neutralizes** bases, most acids and solvents.

How to use it?

Act as fast as possible, preferably less than 10 seconds after the splash

- Open the bottle by turning the lid,
- Gently pour the contents of the bottle into the eye while keeping it open,
- Ask colleagues to fetch for another bottle,
- Call **115** ... A first aid will come with a larger bottle and keep rinsing for 15 minutes,
- If necessary continue rinsing with tap water on the eyeshower from the sink.

Go as soon as possible to the ophthalmic hospital (av de France 15 in Lausanne, tel 021 626 81 11)

Expiry date?

The vials must be changed every two years.

Next change January 2020.

GB déc 2017